

Quick Installation Guide

Dino AM Series MAC installation

Dino-Lite Microscope Mac installation:

1. Insert Driver CD
2. Click Dino > Mac> Your model (eg. AM411T series open AM411T\_AM413T)
3. Click Dino-Lite & Dino-Eye Driver OSX 1.01.0001.dmg
4. Click Dino-Lite & Dino Driver Install
5. When an Install Window appears, click "Continue"
6. When "Import Information" window appears, click "Continue"
7. When "Select a Destination" window appears, click "Continue"
8. When the "Easy Install on Untitled" window appears, click "Install"
9. Key in your password if requested
10. When "Install Software" window appears , click "Close"
11. Now restart your computer
12. Connect the Dino-Lite digital Microscope to USB port
13. Double click "DinoScope" icon on your display
14. Your DinoScope Program should now run.

Dino-Lite support for MAC OS									
Minor version	.0	.1	.2	.3	.4	.5	.6	.7 /.8/.9	.10/.11
X10.0	No	No	No	No	No				
X10.1	Ok	Ok	Ok	Ok	Ok	Ok			
X10.2	Ok	Ok	Ok	Ok	Ok	Ok	Ok	Ok	
Major version									
X10.3	Ok*	Ok*	Ok*	Ok*	Ok*	Ok*	Ok*	Ok*	
X10.4	Ok	Ok	Ok					Ok	Ok
X10.5 (Leopard)	Ok	Ok							
9	No	Ok	Ok						

(\*) QuickTime 6.5.1 or later is required.  
 AM211/AM310 not Support for MAC OS 10.4.7  
 Tiger OS 10.4.10 / .11 - ok  
 Leopard OS 10.5.0/ .1 - ok